



ULTRA LOW FREEZER MV85 SERIES

INSTRUCTIONS

USER MANUAL

WARNING: READ BEFORE CONTINUING

To reduce the risk of fire, electric shock or injury to persons using this freezer, read all instructions and follow basic safety precautions before using the unit, including the following:



Do not modify the plug provided with the freezer. If it will not fit the outlet, have a proper outlet installed by a qualified electrician.



Do not position equipment so it is difficult to disconnect from the power supply. freezer must be at least 6" away from any wall or object on any side.



While under warranty, do not attempt to repair or replace any part of the freezer for servicing without first contacting the So-Low Service Department.

SAVE THESE INSTRUCTIONS

So-Low Environmental Equipment Company
10310 Spartan Drive
Cincinnati, OH 45215-1221
Tel: 513-772-9410
<http://www.so-low.com>

For customer service:
Email: sales@so-low.com

For parts replacement:
Email: parts@so-low.com

For technical support:
Email: service@so-low.com

CONTENTS

◆ SYMBOLS AND STARTING INSTRUCTIONS.....	1
• MEANING OF ILLUSTRATED SYMBOLS.....	1
• STARTING INSTRUCTIONS.....	1
◆ UNIT REQUIREMENTS.....	2
• PRE-INSTALLATION INFORMATION.....	2
◆ GENERAL SPECIFICATIONS.....	3
• TEMPERATURE SPECIFICATIONS.....	3
• ELECTRICAL SPECIFICATIONS.....	3
◆ REFRIGERATION SPECIFICATIONS.....	4
◆ MAINTENANCE.....	5
◆ UNIT OPERATION.....	6
◆ TEMPERATURE CONTROL.....	7
◆ ALARM OPERATION.....	8
• ALARM SYSTEM.....	8
◆ CONTROL PROGRAMMING.....	9
• CUSTOM SETTINGS.....	9
• ADVANCED SETTINGS.....	10
• SERVICE SETTINGS.....	11
• ADVANCD SERVICE SETTINGS.....	11
◆ DATA DOWNLOAD.....	12
• DOWNLOAD.....	12
• UPLOAD.....	12
◆ FACTORY SETTINGS.....	13
◆ WIRING DIAGRAM.....	14
• MV85-1 DIAGRAM.....	15
• MV85-1 PARTS LIST.....	16
• MV85-2 DIAGRAM.....	17
• MV85-2 PARTS LIST.....	18
• MV85-3 DIAGRAM.....	19
• MV85-3 PARTS LIST.....	20

SYMBOLS AND STARTING INSTRUCTIONS

Explanation

MEANING OF ILLUSTRATED SYMBOLS

ILLUSTRATED SYMBOLS

Various symbols are used in this safety manual in order to use the unit without danger of injury and damage of the unit. A list of problems caused by ignoring the warnings and improper handling is divided as shown below. Be sure that you understand the warnings and cautions in this manual before operating the unit.



CAUTION

BLACK WITH YELLOW BACKGROUND
LIGHTNING BOLT
CAUTION, RISK OF ELECTRICAL SHOCK



WARNING

BLACK WITH YELLOW BACKGROUND
EXCLAMATION POINT
CAUTION, REFER TO ACCOMPANYING DOCUMENTS

STARTING INSTRUCTIONS

1. Move the freezer to an indoor location, and plug the freezer into an appropriate outlet with an adequate power supply. Consult your maintenance department for additional information on the proper electrical configuration for this unit.
2. Once plugged in, the compressor(s) will start to operate and pull down to the setpoint on the temperature control.
3. Allow the freezer to reach the setpoint temperature. Depending on the size of the unit, this may take up to 12 hours.
4. Product can now be loaded into the freezer for storage.

Unit Information

PRE-INSTALLATION INFORMATION

RANGE OF ENVIRONMENTAL CONDITIONS FOR WHICH THIS EQUIPMENT IS DESIGNED

1. Indoor use.
2. Altitude up to 2000m.
3. Ambient temperatures 15°C to 30°C (60°F TO 85°F)
4. Recommended humidity range of 30% to 90%.
5. Mains supply fluctuations up to -5% to +10% of the nominal voltage.
6. Transient over-voltages typically present on the mains supply (overvoltage category II). Pollution degree 1.



WARNING

THIS FREEZER IS PROVIDED WITH AN INPUT CIRCUIT PROTECTIVE DEVICE WHICH SHALL BE MAINTAINED AND SERVICED BY QUALIFIED PERSONNEL ONLY.



CAUTION

UNDER NO CIRCUMSTANCES USE EXTENSION LEADS OR ADAPTORS. CONNECT THE APPLIANCE DIRECTLY TO A FIXED INSTALLATION. IF THE APPLIANCE IS CONNECTED TO A LONG EXTENSION CABLE THERE IS A RISK THAT THE CABLE WILL BECOME HOT.



CAUTION

UNPLUG FREEZER BEFORE ANY TECHNICAL SERVICE IS PERFORMED ON THE UNIT.



WARNING

DO NOT POSITION EQUIPMENT SO IT IS DIFFICULT TO DISCONNECT FROM THE POWER SUPPLY.

GENERAL SPECIFICATIONS

Specifications

TEMPERATURE SPECIFICATIONS


OPERATIONAL TEMPERATURE RANGE

-40°C TO -85°C

ELECTRICAL SPECIFICATIONS

ELECTRICAL PLUG

- Plug the freezer into the proper outlet with an adequate power supply.
- This unit requires a **Dedicated Electrical Line**.

MODEL	VOLTAGE	AMPERAGE	PLUG
MV85-1 MV85-2 MV85-3	115 VOLTS 60 HERTZ 1 PHASE	15 AMP DEDICATED LINE	 NEMA 5-15



WARNING

ONLY PLUG THIS UNIT INTO THE PROPER OUTLET. DO NOT ATTEMPT TO MODIFY PLUG IN ANY WAY. IMPROPER USE OF THE ELECTRICAL PLUG WILL VOID WARRANTY



CAUTION

FAILURE TO GROUND THE EQUIPMENT MAY CAUSE PERSONAL INJURY OR DAMAGE TO THE EQUIPMENT. ALWAYS CONFORM TO THE ELECTRICAL CODES.

REFRIGERATION SPECIFICATIONS

REFRIGERATION SPECIFICATIONS

MODEL	CU. FT.	REFRIGERATION SYSTEM	REFRIGERANT
MV85-1	0.25	3/8 HP REFRIGERATION SYSTEM	EP88 – 6.10 OZ.
MV85-2	2	3/8 HP REFRIGERATION SYSTEM	EP88 – 6.88 OZ.
MV85-3	3	3/8 HP REFRIGERATION SYSTEM	EP88 – 6.90 OZ.

PREVENTATIVE MAINTENANCE

BEFORE PERFORMING MAINTENANCE

To reduce the risk of fire, electric shock or injury to persons using this freezer, read all instructions and follow basic safety precautions.



CAUTION

DISCONNECT THIS UNIT FROM THE POWER SUPPLY BEFORE PERFORMING MAINTENANCE ON THE UNIT.

CLEANING PROCEDURE

- Wipe down the exterior of the freezer with a soft cloth and spray type polish.
- If frost builds up in the chamber, a bucket and ice-scraper can be used to remove the ice.
- If excessive ice builds up, the unit can be defrosted (see below).

DEFROST PROCEDURE

1. Remove any product in the freezer and store it in a back-up freezer or elsewhere.
 2. Unplug the freezer, and open the freezer front door / lid.
 3. Use a cloth to protect the control from dripping water.
 4. Air out the freezer for at least 12 hours, allowing the unit to reach room temperature.
 5. Take a rag and wipe up all the excess water in the unit (melted frost).
 6. Plug the unit in and set your temperature to the desired setpoint
 7. Once the desired temperature is reached, add product back into the unit.
-
-



WARNING

IT IS RECOMMENDED TO SLOWLY RE-ADD YOUR PRODUCT INTO THE FREEZER TO PREVENT AN EXTREME LOAD ON THE COMPRESSORS, WHICH COULD SHORTEN FREEZER LIFE EXPECTANCY.

GENERAL OPERATION

- The freezer should be located in a dry, cool place without direct sunlight.
- To start the freezer, set the digital thermometer at the temperature required. It will now take about 6-10 hours before the set temperature is reached.
- The freezer should be cooled down empty.



WARNING

WHEN FILLING THE UNIT WITH RACKS/BOXES, MAKE SURE TO PRE-COOL THEM IN ORDER NOT TO EFFECT THE INSIDE TEMPERATURE OF THE UNIT.



WARNING

THE INNER DOORS SHOULD BE USED AT ALL TIMES. MAKE SURE TO CLOSE THEM TIGHTLY.

THE MANUFACTURER IS NOT RESPONSIBLE FOR DAMAGE TO THE FREEZER CAUSED BY FAILURE TO DO SO.




WARNING

CLEAN THE FREEZER BOTH INSIDE AND OUT USING A DAMP, WRUNG OUT CLOTH. DRY WITH A DRY CLOTH.

THE SMELL FROM THE PLASTIC PARTS IN THE FREEZER WILL DISAPPEAR ONCE THE FREEZER HAS BEEN COOLED DOWN.


G-214 CONTROLLER

CONTROL DISPLAY

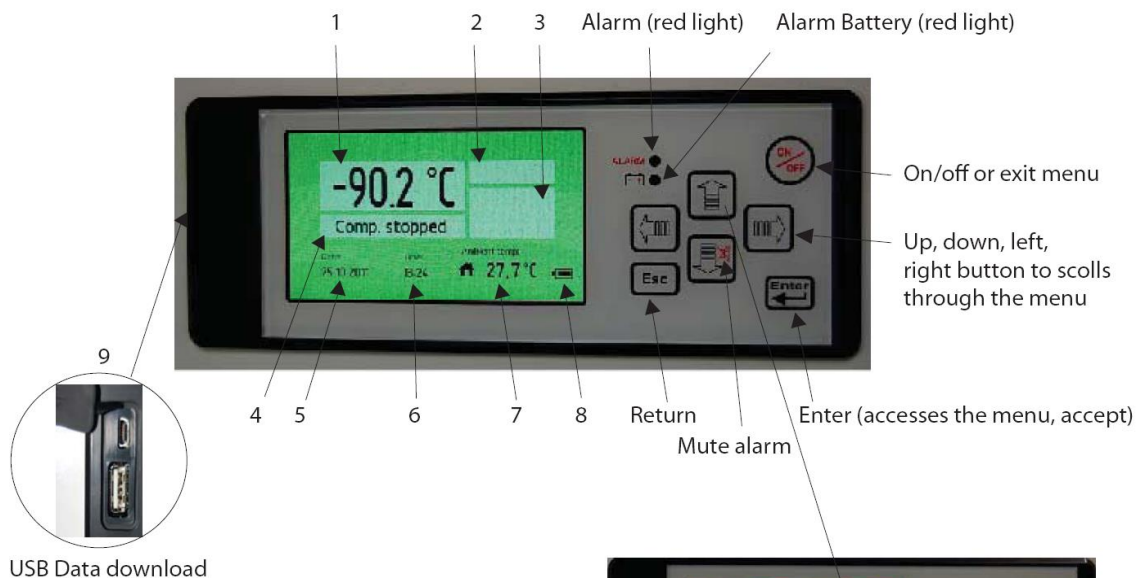

Control Panel Description

1. Temperature indicator
2. Alarm indicator
3. Alarm-icons
4. Compressor indication
5. Date indication
6. Hour indication
7. Ambient temperature indicator
8. Battery level
9. USB-Connection

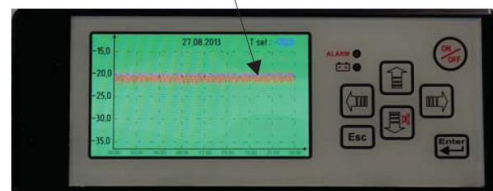

Push the **UP** button in order to generate a graph.

Push the **LEFT** button in order to return up to 10 days back.

Push the **Esc** button in order to exit the menu.

ALARM SYSTEM

Alarms on display



Alarm icons



Door open alarm

The **Door open alarm** indicate, that the door is not correct closed



High temp. alarm

The **High temperature alarm** indicate, that the temperature inside the unit is higher than permitted from setting of the unit.



Low temp. alarm

The **Low temperature alarm** indicate, that the temperature inside the unit is lower than permitted from setting of the unit.



Probe eprom failure

The **Probe eprom failure** alarm indicate, that the probe is not working.



Power failure


The **Power failure** alarm indicate, that there is no power to the unit.




Low battery

The **Low battery** alarm indicate, that the power on the battery is to low.

CUSTOM SETTINGS

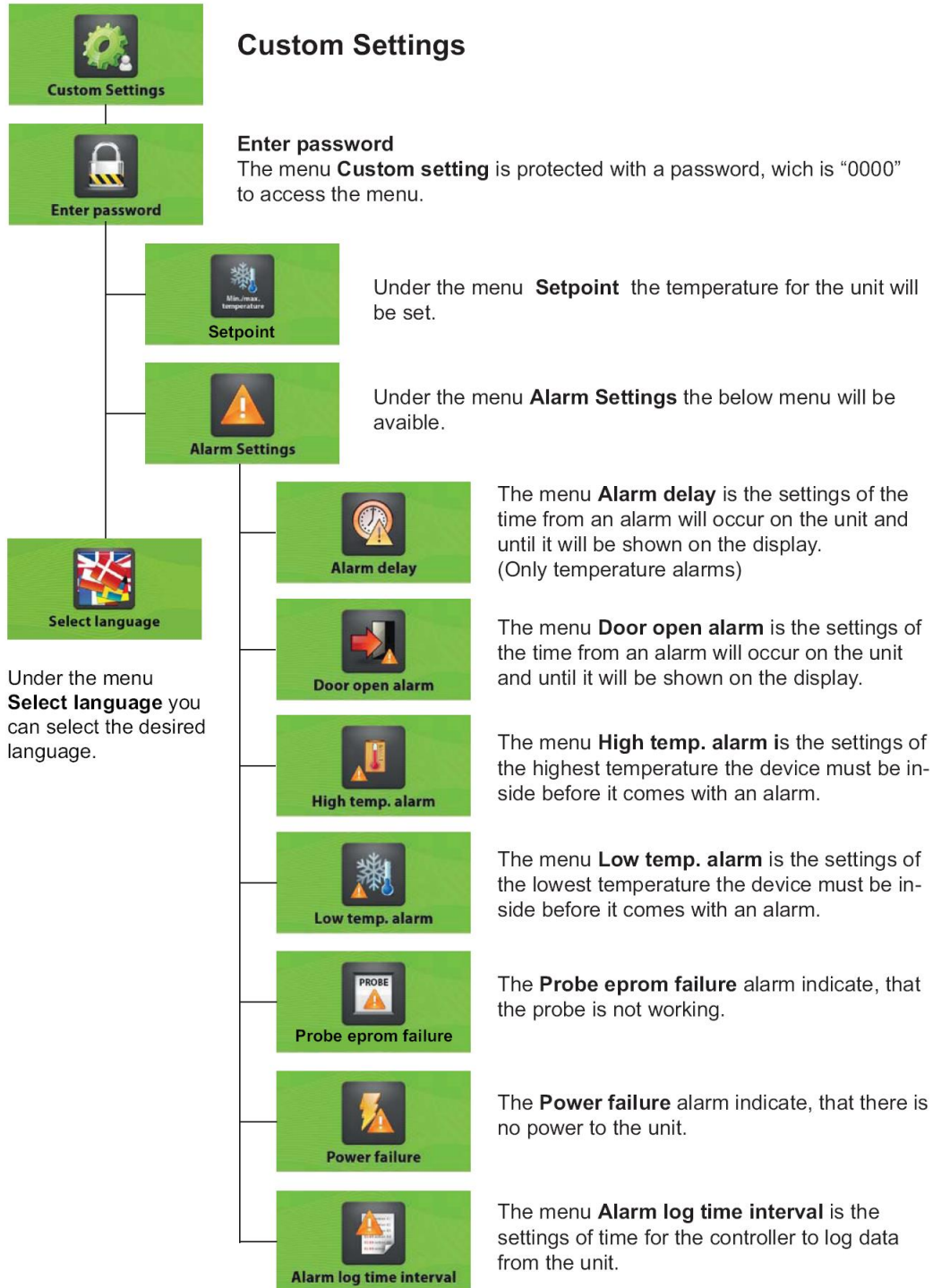

ADVANCED SETTINGS

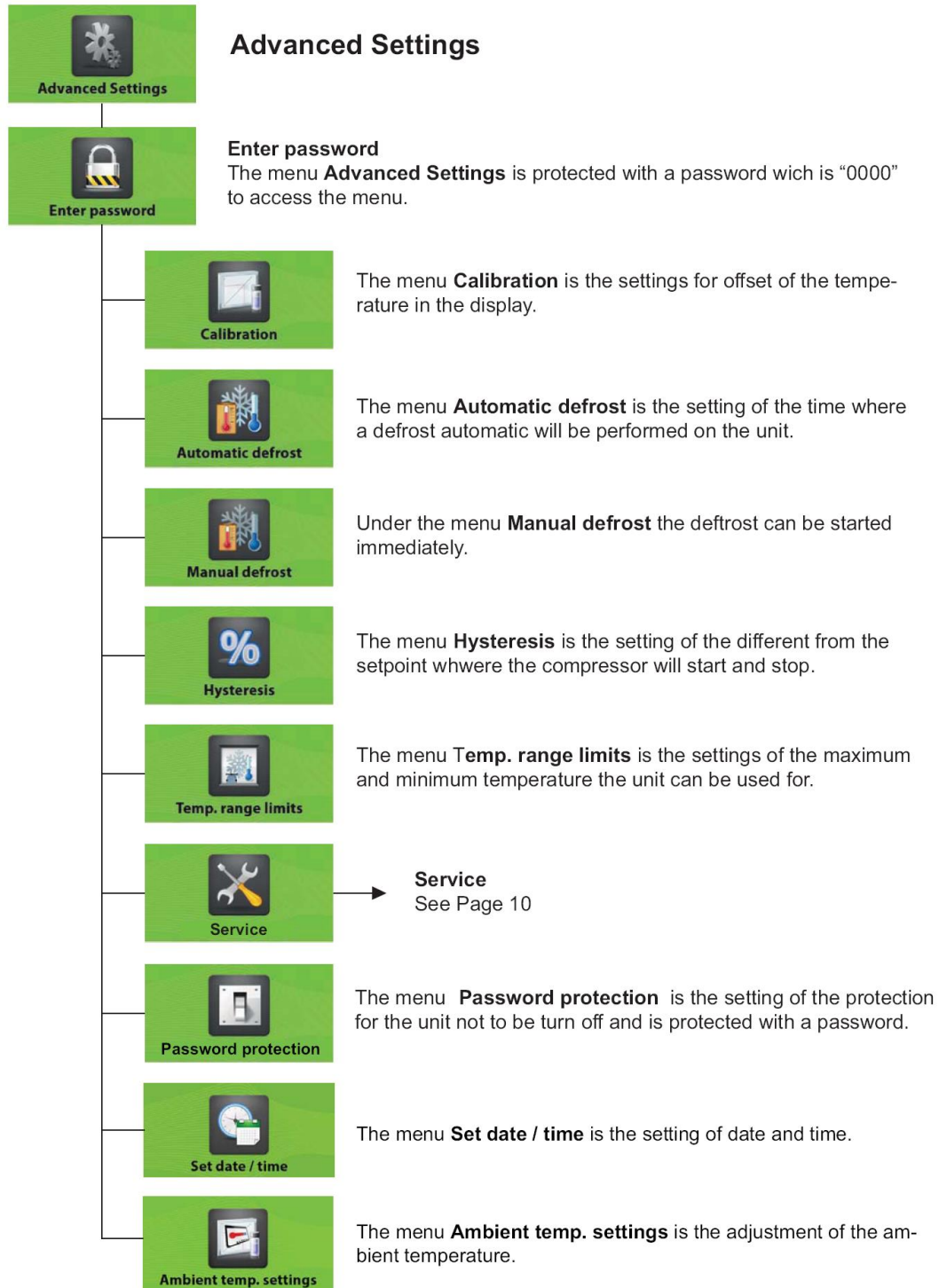

SERVICE SETTINGS



Service

Service Settings

Under the menu **Service**, there is information about the unit, which is important for service on the unit.



Compressor hours

The menu **Compressor hours** shows, how many hours the compressor has run.



Fan hours

The menu **Fan hours** shows, how many hours the fan has run.



Probe

Under the menu **Probe** the temperature for the sensor mounted in the unit is displayed.



Probe inside unit

Under the menu **Probe inside unit** the temperature of the sensor mounted inside the unit is displayed. There is mounted a sensor as standard in the unit, but it is possible to mount 3 sensors inside the device. (Controller version)



Probe Compressor

Under the item **Probe Compressor** the temperature of the sensor mounted on the compressor is displayed. (Not all models)



Probe evaporator

Under the point **Probe evaporator** the temperature of the sensor, that is located on the condenser, is displayed.

ADVANCED SERVICE SETTINGS



Advanced Service Settings

Advanced Service Settings

This menu is only used by the service department at the manufacturer.

DOWNLOAD


Download Data



Place the USB pen drive in the connection on the display and press the button right for **read data** from the unit to the USB pen.



The **Read data** will be shown in the display when the data is loading to the USB pen drive.

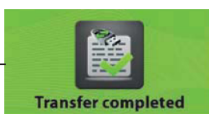

The **Transfer complete** will be shown on the display when the loading is finished and the USB pen can be removed from the display.


Open the 2 files:
data00 and **param00**
on a computer in Excel or similar.

UPLOAD


Upload Data (Only for new settings from the manufactory)



Turn-off the display before start.
Place the USB pen drive in the connection on the display and press the button left for upload parameters from the USB pen to the unit.

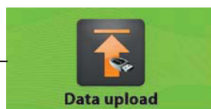

The **Data upload** will be shown in the display when the data is loading to the unit.

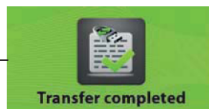

The **Transfer complete** will be shown on the display when the loading is finished and the USB pen can be removed from the display.

FACTORY SETTINGS


BASIC SETTINGS

DESCRIPTION	DEFAULT SETTINGS	CUSTOMER SETTING
PASSWORD	0000	
TEMPERATURE SETPOINT	-80.0	
ALARM – DELAY	15 Min	
ALARM – DOOR OPEN	Enable	
ALARM – HIGH TEMPERATURE	-70	
ALARM – LOW TEMPERATURE	-90	
ALARM – LOG TIME INTERVAL	1	
PROBE (EPROM) FAILURE	Enable	
POWER FAILURE	Enable	


ADVANCED SETTINGS

DESCRIPTION	DEFAULT SETTINGS	CUSTOMER SETTING
PASSWORD	0000	
CALIBRATION	0.0	
AUTOMATIC DEFROST	0	
MANUAL DEFROST	OK	
HYSTERESIS	2.0	
MAXIMUM RANGE LIMIT	-40	
MINIMUM RANGE LIMIT	-86	

WIRING DIAGRAM

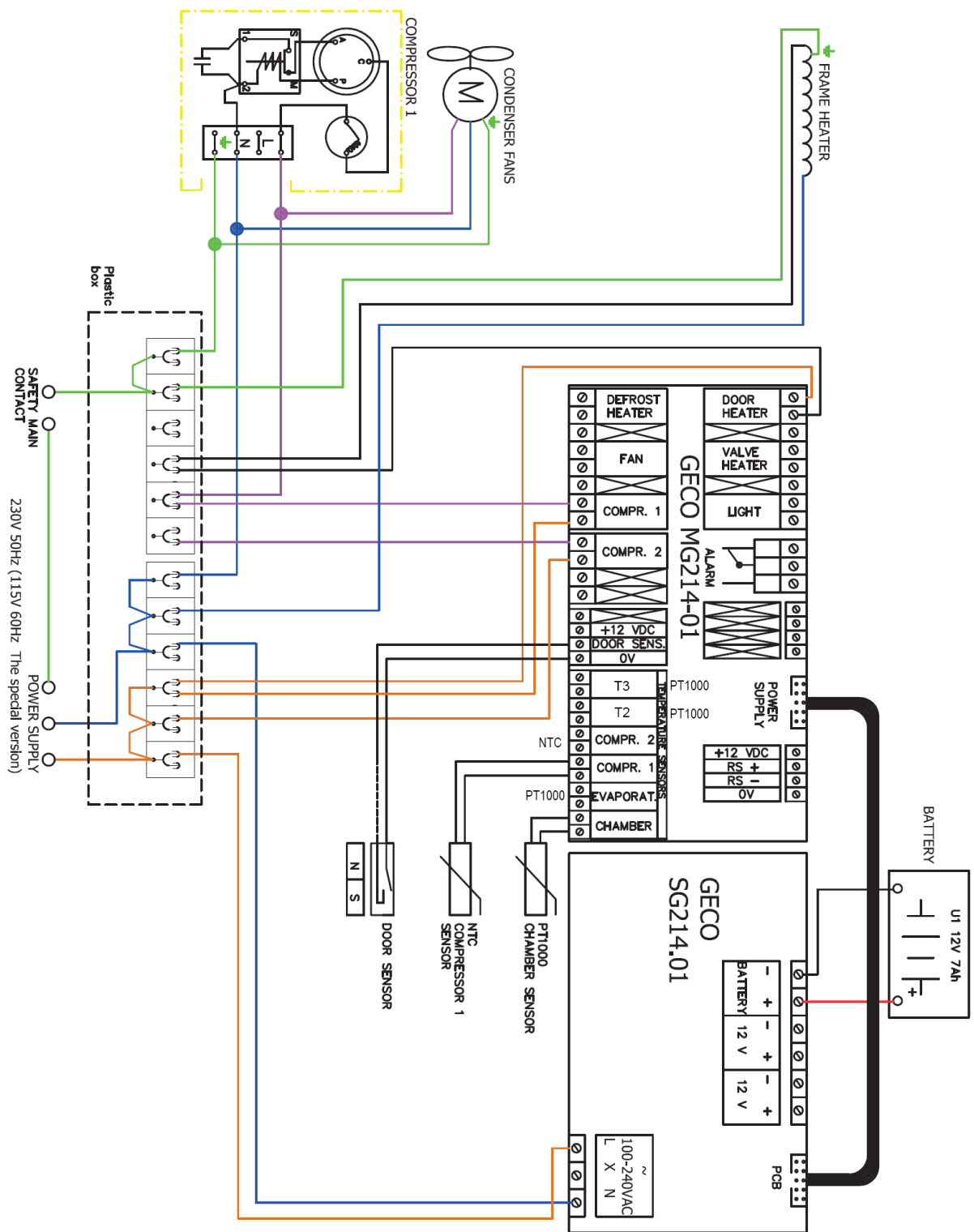


MV85-1 DIAGRAM

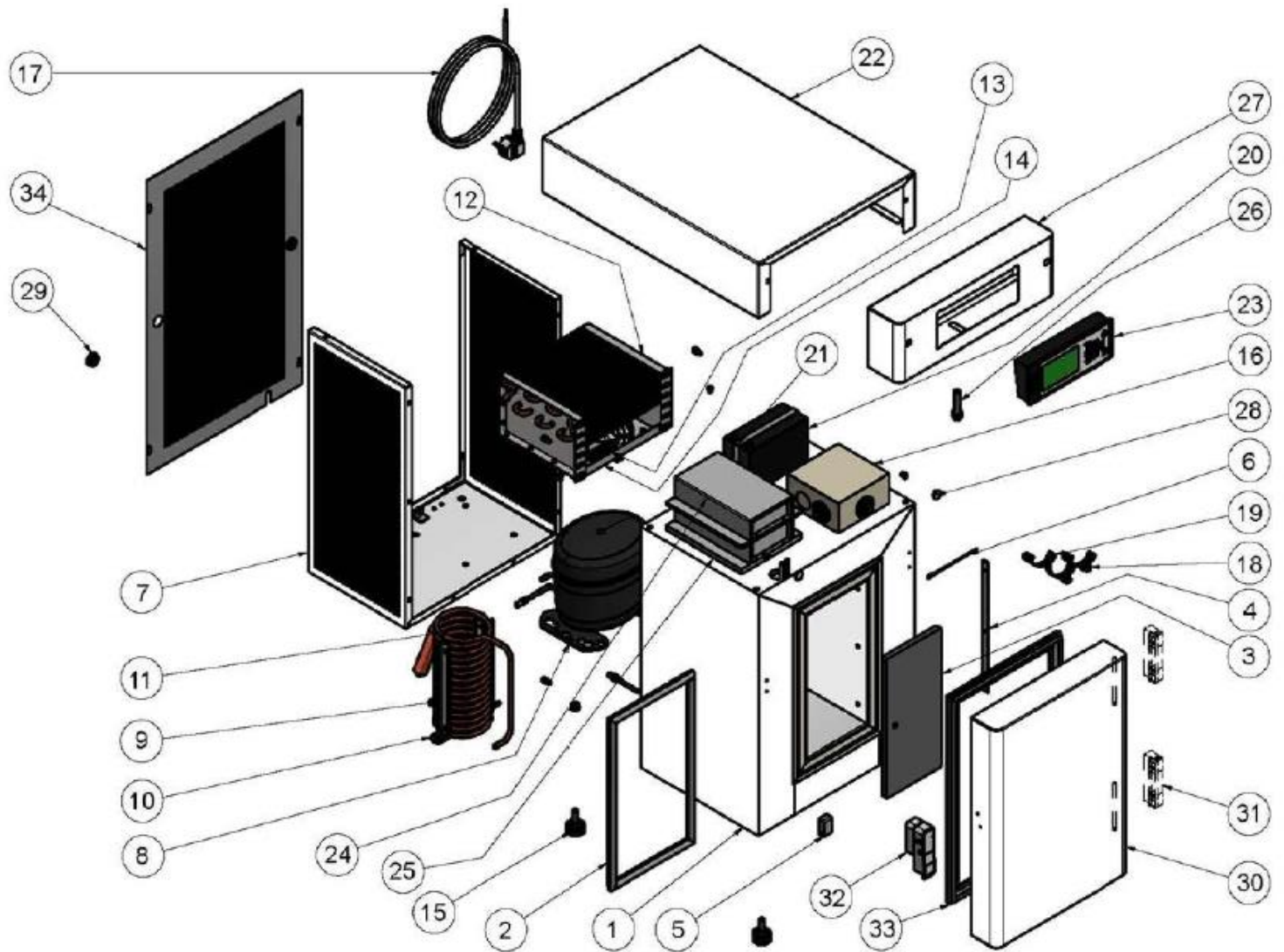

MV85-1 PARTS LIST

Pos. no.	Code no.	Part	Recommended
1	-	Cabinet	
2	5052159-01	Gasket for frame (silicone)	
3	5052136-01	Inner door	
4	5031218-01	Hinge for inner door - 280mm	
5	5051065-01	Handle (Belt) for inner door	
6	5042003-01	Probe (PT1000)	x
7	5031228-01	Aggregate cover	
8	5070007-01	Compresseor SC12CNX.2	
9	5074003-01	Condenser spiral	
10	5031234-01	Condenser spiral support 1	
11	5031235-01	Condenser spiral support 2	
12	5071039-01	Condenser TK0680	
13	5010007-01	Condenser fan ex.wing	x
14	5010043-01	Fan wing	
15	5053030-01	Regulated feet	
16	5049006-01	Electrical box for 230V connection G-214	
17	5019169-01	Mains lead ULUF	
18	5019128-01	Flat cable G-214 PCB and PSU	
19	5019017-01	Flat cable G-214 display	
20	5044001-01	Battery G-210/214 controller	x
21	5031192-01	Battery support	
22	5031230-01	Top cover ULUF 15	
23	5043007-01	Display for G-214	x
24	5041007-01	Power supply G-214	x
25	5041006-01	PCB-board G-214	x
26	5019113-01	Door switch	
27	5031231-01	Front panel	
28	5030098-01	Hole plug for cabinet	
29	5030099-01	Hole cover for back cover	
30	5052156-01	Door ULUF 15	
31	5031025-01	Hinges	
32	5051006-01	Lock A-458-R	
33	5052158-01	Door magnet. Gasket ULUF 15	
34	5031229-01	Back cover	
-	5079000-01	Service dryfilter	x
-	5072000-01	Cylinder + deposit	x
-	5072023-01	Refrigerant EP88 - 173g	x

MV85-2 DIAGRAM

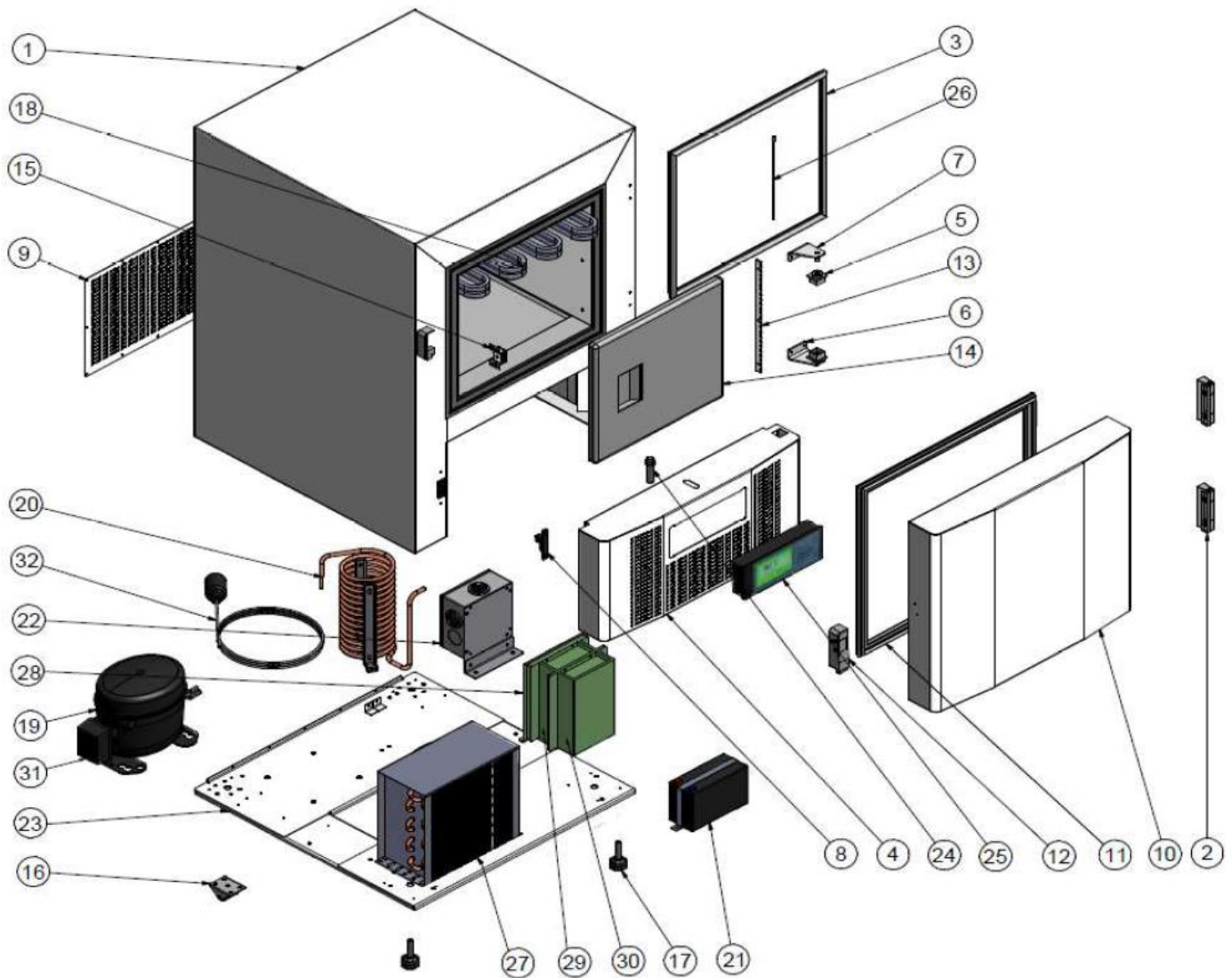

MV85-2 PARTS LIST

[illegible]

This exploded view diagram illustrates the assembly of a microwave oven. The components are numbered as follows:

- 1: Microwave cabinet body
- 2: Top panel
- 3: Control panel
- 4: Control panel mounting bracket
- 5: Control panel mounting screw
- 6: Control panel mounting screw
- 7: Control panel mounting screw
- 8: Control panel mounting bracket
- 9: Control panel mounting screw
- 10: Control panel mounting screw
- 11: Microwave door
- 12: Door hinge
- 13: Door hinge
- 14: Door hinge
- 15: Door hinge
- 16: Door hinge
- 17: Door hinge
- 18: Door hinge
- 19: Door hinge
- 20: Door hinge
- 21: Door hinge
- 22: Door hinge
- 23: Door hinge
- 24: Door hinge
- 25: Door hinge
- 26: Door hinge
- 27: Door hinge
- 28: Door hinge
- 29: Door hinge
- 30: Door hinge
- 31: Door hinge
- 32: Door hinge

MV85-3 PARTS LIST

Pos. no.	Code no.	Part	Recommended
1		Cabinet	
2	5031153-01	Side cover	
3	5050053-01	Front panel	
4	5031157-01	Back panel hinge	
5	5031154-01	Back panel	
6	5030082-01	Hole cover ø22	
7	5031025-01	Set of hinges	
8	5052123-01	Door, complete	
9	5052125-01	Door magnetic gasket	
10	5051006-01	Latch incl. Set of keys	
11	5052127-01	Silicone gasket	
12	5049005-01	Door switch	
13	5031150-01	Subdoor hinge	
14	5060024-01	Shelf	
15	5052129-01	Subdoor	
16	5031013-01	Magnetic latch for subdoor	
17	5073033-01	Evaporator	
18	5031151-01	Electrical support	
19	5044001-01	Battery BP 7-12 (12V, 7AH)	1
20	5031139-01	Power supply support	
21	5049006-01	Electrical box for 230V connection	
22	5071046-01	Condenser TK1000	
23	5070015-01	Compressor SC18CNX.2 Complete	
24	5074003-01	Condenser spiral	
25	5043007-02	Display G-214	1
26	5053029-01	Castor	
27	5053016-01	Regulated feet M10x40	
28	5031156-01	Agregate base	
29	5041007-01	Power supply G-214	1
30	5041006-01	PCB Board G-214 Controller	1
31a	5070021-01	Starting capacitor 117U5017	1
31b	5070026-01	Starting device 117U7011	1
32	5042003-01	Probe (PT 1000)	1
33	5019169-01	Main Lead	
-	5042016-01	Probe NTC for Compressor	1
-	Finans no.	Refrigerant 195g EP88	1
-	5072000-01	Cylinder deposit for refrigerant	1
-	5079000-01	Service dryfilter	1